

Diaphragm monitoring system with HART[®] protocol For sanitary applications Model DMSU21SA



for further approvals see page 9

Applications

- Hygienic pressure measurement for the pharmaceutical industry and for aseptic food processing
- Pressure/vacuum measurement on pipelines, fermenters, bioreactors and vessels as well as with processing and transport of high-quality media
- Suitable for the production of active pharmaceutical ingredients (API)
- For the monitoring of processes with sterile steam
- For gases, vapour; liquid, pasty, powdery and crystallising media

Special features

- Double-diaphragm system prevents contamination of process and environment
- Hygienic process connections in different designs
- Signal transmission and configuration with only one cable per measuring location
- Minimal installation costs, even with retrofitting

Description

The model DMSU21SA diaphragm monitoring system fulfils the constantly growing requirements for process monitoring on plants for sanitary applications. With the HART[®] protocol, not only the measuring signal, but also the status of the integrated diaphragm monitoring can be transmitted, as required, to the process control. The system architecture with integrated diaphragm monitoring bridges the gap between digitalisation and process reliability.

With the patented diaphragm monitoring, a monitoring element is mainly used for electrical/digital signal transmission of the diaphragm condition. In addition, the diaphragm condition is displayed on a dial with red/green areas. In the event of a diaphragm rupture, a second, internal diaphragm ensures the reliable separation of the environment and the process. A fluid inside the system, which is explicitly chosen to suit the particular measuring requirement, hydraulically transmits the pressure to the pressure measuring instrument. Depending on the application, the system fill fluid is FDA- and also USP-compliant.

The diaphragm monitoring system can be supplied with all common process connections for sanitary applications and fulfils the applicable standards for them, e.g. EHEDG, 3-A and ASME BPE.

This product, WIKA model DMSU21SA, is a patented WIKA development (US 10794787, NL 2019251).

WIKA data sheet DS 95.11





Diaphragm monitoring system, model DMSU21SA

WIKA data sheet DS 95.11 · 05/2021


Data sheets showing similar products:

Diaphragm monitoring system with clamp connection; model DMS-FP; see data sheet DS 95.20 Diaphragm monitoring system with threaded connection; model DMS34; see data sheet DS 95.18 Diaphragm pressure gauge, flush; with integrated diaphragm element monitoring; model PG43SA-D; see data sheet PM 04.17 Process transmitter, with pressure port or flush diaphragm; models UPT-20, UPT-21; see data sheet PE 86.05 Page 1 of 27




Illustrative example of a diaphragm monitoring system

The diaphragm monitoring system is a combination of pressure measuring instrument and diaphragm seal with an additional monitoring element for the diaphragm condition.

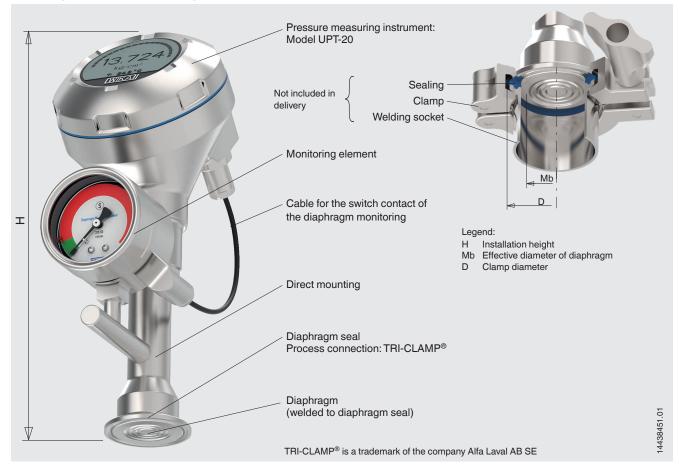

Functional principle of diaphragm monitoring

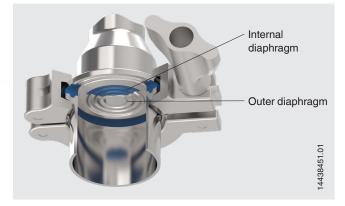

Diaphragm rupture

In the event of a diaphragm rupture, the pressure monitored in the intermediate space increases. As soon as the display of the monitoring element exceeds the predefined set point, the alarm signal for a diaphragm rupture is transmitted. The alarm signal is output either via the HART[®] protocol or as an error signal on the current loop. The simple cabling enables a quick instrument replacement at the measuring location. WIKA's double-diaphragm design is the solution for critical processes where neither the medium should find its way into the environment, nor should the system fill fluid find its way into the product.

Normal operation

In normal operation, the pressure measurement and the diaphragm monitoring operate without limitations within the performance limits of the overall system.

The space between the two diaphragms is evacuated. With the monitoring element, this vacuum is measured and the condition signalled in the green area; no electrical/digital alarm signal will be output.

Safety

The measuring technology of the monitoring element withstands the process pressure despite the diaphragm rupture. The measuring function of the overall system is maintained without limitations. The process safety is guaranteed because the materials used for the two diaphragms are the same as those of the wetted parts of the diaphragm seal. The overall system is, nevertheless, damaged and must be replaced immediately.

Specifications

Diaphragm seal system	
Design	Model UPT-20 process transmitter mounted to a diaphragm seal with clamp connection, welded
Material ¹⁾	
Wetted	Diaphragm and diaphragm seal: Stainless steel 1.4435 (316L); UNS S31603
Surface roughness	
Wetted	 Ra ≤ 0.38 μm [15 μin] per ASME BPE SF4, electropolished (except for weld seam) Ra ≤ 0.76 μm [30 μin] (except for weld seam)
Non-wetted	$Ra \le 0.76 \ \mu m [30 \ \mu in]$ (except for weld seam)
Case head	 Plastic (PBT) with conductive surface per EN 60079-0:2012, colour: Night blue RAL5022 Stainless steel case 1.4308 (CF-8), precision-cast (suitable for chemical and petrochemical industries) Stainless steel case 1.4308 (CF-8) with electropolished surface (suitable for pharmaceutical, food and hygienic industries)
System fill fluid	 Medicinal white mineral oil, FDA 21 CFR 172.878 Medicinal white mineral oil, FDA 21 CFR 178.362 (a)
	USP-, EP- and JP-certified USP = United States Pharmacopeia EP = European Pharmacopoeia JP = Japanese Pharmacopoeia
Mounting type	Direct mounting
Level of cleanliness of wetted parts	Oil- and grease-free per ASTM G93-03 level F WIKA standard (< 1,000 mg/m ²)

1) Other materials on request

Diaphragm monitoring ¹⁾ via switch contact and dial with red/green areas			
Output signal	The status of the switch contact (alarm signal) is monitored in the process transmitter and output via the HART [®] protocol or as an error signal on the current loop. \rightarrow see "Output signals"		
Alarm signal condition	For the monitoring element to change the status of the switch contact, the specified set point must be reached for at least 1.5 seconds. This prevents shocks or vibrations from triggering the alarm signal unintentionally.		
Dial display	Pointer in green range \rightarrow Outer diaphragm intactPointer in red range \rightarrow Outer diaphragm defective		
Material			
Case	Stainless steel, with solid baffle wall (Solidfront) and blow-out back		
Pressure element	Stainless steel 316L		
Movement	Stainless steel		
Bayonet ring	Stainless steel		
Pointer/dial	Aluminium		
Window	Laminated safety glass		

1) In the event of a diaphragm rupture, the diaphragm monitoring is specified for process pressures ≥ 0 bar gauge

Measuring range

Gaug	e pressure					
bar	0 2.5	04	06	0 10	0 16	0 25
psi	0 30	0 60	0 100	0 160	0 250	0 360

Vacuum and +/- measuring ranges ¹⁾						
bar	-1 +1.5	-1 +3	-1 +5	-1 +9	-1 +15	-1 +24
psi	-14.5 20	-14.5 +40	-14.5 +80	-14.5 +130	-14.5 +200	-14.5 +350

1) Under vacuum, diaphragm monitoring is only possible to a limited extent

Other units settable (\rightarrow see page 7).

Other measuring ranges can be set via turndown.

For example, a 0 ... 6 bar [0 ... 100 psi] instrument can also be used from -1 ... +6 bar [-14.5 ... +100 psi].

Vacuum/overload safety		
Vacuum safety	Yes	
Overload safety	1-fold	

Output signal

Output signal	
Signal types	 4 20 mA with HART[®] signal (HART[®] rev. 7) 4 20 mA
Load in Ω	\leq U+ - 14 V / 0.023 A
	U+ = Applied supply voltage (→ see "Supply voltage")
Dampening	0 99.9 s, adjustable
	After the set dampening time the instrument outputs 63 % of the applied pressure as output signal.
Settling time t ₉₀	80 ms
Refresh rate	50 ms

Process connection

Standard	
Clamp connection per DIN 32676	→ See from page 11
Clamp connection per ISO 2852	→ See page 13
Aseptic threaded pipe connection per DIN 11864-1	→ See from page 14
Aseptic flange connection per DIN 11864-2	→ See from page 17
Aseptic clamp connection per DIN 11864-3	→ See from page 20
Ingold connection with union nut	→ See page 23

Accuracy specifications

Accuracy specifications			
Accuracy at reference conditions	 0.1 % of span 0.5 % of span 		
Adjustability			
Zero point	-20 \ldots +95 % (downwards, the adjustability is always limited by the minimum pressure of 0 bar abs. [0 psia])		
Span	-120 +120 $\%$ with a difference between zero point and span of max. 120 $\%$ of the nominal measuring range		
Turndown	Unlimited; maximum recommended turndown 20:1		
	Measuring range ≤ 25 bar [360 psi]		
Mounting correction	-20 +20 %		
Non-repeatability	\leq 0.1 % of span		
Behaviour with turndown ²⁾			
TD ≤ 5:1	No influence on the accuracy		
TD > 5:1 ≤ 100:1	GES = GG x TD / 5		
Long-term stability	\leq 0.1 % of span		

1) Including non-linearity, hysteresis, zero offset and end value deviation (corresponds to measured error per IEC 61298-2).

2) Legend GES: Overall accuracy via turndown GG: Accuracy (e.g. 0.15 %) TD: Turndown factor (e.g. 4:1 corresponds to TD factor 4)

Reference conditions per IEC 61298-1

Reference conditions per IEC 61298-1		
Temperature	23 °C ±2 °C [73 °F ±7 °F]	
Supply voltage	DC 23 25 V	
Air pressure	860 1,060 mbar [86 106 kPa, 12.5 15.4 psi]	
Air humidity	45 75 % r. h.	
Characteristic curve determination	Terminal method per IEC 61298-2	
Curve characteristics	Linear	
Reference mounting position	Vertical, diaphragm points downward	

Voltage supply

Voltage supply	
Supply voltage U+	DC 14 30 V
Maximum voltage U _i	DC 30 V
Maximum current li	100 mA
Maximum power P _i (gas)	1,000 mW
Effective internal capacitance	11 nF
Effective internal inductance	100 μH

→ For more information, see "Approvals"

Case design of process transmitter

Case	
Material	 Plastic (PBT) with conductive surface per EN 60079-0:2012. Colour: Night blue RAL5022 Stainless steel case 1.4308 (CF-8) with electropolished surface (suitable for pharmaceutical, food and hygienic industries)


Electrical connection

Electrical connection		
Cable glands M20 x 1.5		
Stainless steel cable gland in hygienic design	Sealing	FDA-compliant
	Cable diameter	6 12 mm [0.24 0.47 in]
	Ingress protection per IEC/EN 60529	IP66/67
Angular connector DIN 175301- 803A with mating connector	Wire cross-section	Max. 1.5 mm ² (AWG 16)
	Ingress protection per IEC/EN 60529	IP65 ¹⁾
Circular connector M12 x 1 (4-pin) without mating connector	Ingress protection per IEC/EN 60529	IP65 ¹⁾
Electrical safety	Reverse polarity protection	


1) The stated ingress protection only applies when plugged in using mating connectors that have the appropriate ingress protection.

Pin assignment

Cable gland M20 x 1.5 and spring-loaded terminals

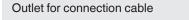

Pin assignment

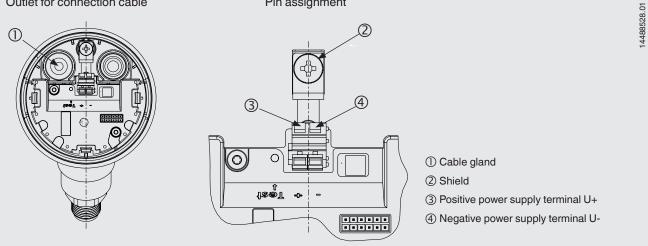

Display and operating unit, model DI-PT-U

Display and operating unit, n	nodel DI-PT-U				
Display type	LC display ¹⁾				
Refresh rate	200 ms				
Main display	4 ½-digit				
Segment display	Character size 14 mm [0.55 in]				
Additional display	Selectable via menu, three-line scale range				
Bar graph display	20 segments, radial, pressure gauge simulation				
Adjustable units					
Pressure units	barftH2OpsiinH2OmbarmHgMPammHgkPainHghPakg/cm²Pag/cm²mH2OTorrmmH2O				
Level units	 m cm mm ft in 				
Volume units	 I m³ gal inch³ ft³ % 				
Free unit	A free character string (6 characters) can be defined as a unit				
Operating state	Display via symbols				
Operating temperature	By using the display and operating unit, the operating temperature is limited to -20 +60 $^\circ C$ [-4 +140 $^\circ F$]				
Storage temperature	-40 +80 °C [-40 +176 °F]				
Operation	4 buttons for the recall and operation of the settings				
Colours					
Background	Light grey				
Digits	Black				
Display size	Separated into main and segment displays				
Ingress protection per IEC/EN 60529	 IP20 (loose) IP40 (built-in without cover) 				
Material	Case from ABS, window from polyester film				

1) For the process transmitter only this display may be used. \rightarrow For order number see "Accessories and spare parts".

Operating conditions

Operating conditions					
Range of applications	Suitable for internal and external operation, exposure to direct sunlight is permitted				
Permissible air humidity	≤ 93 % r. h.				
Permissible temperature ranges					
Ambient	10 40 °C [50 104 °F]				
Medium	 -10 +130 °C [14 266 °F] -10 +150 °C [14 302 °F] 				
Storage	10 60 °C [50 140 °F]				
Ingress protection per IEC/EN 60529	IP65				
	The ingress protection only applies with closed case head and closed cable glands.				
Explosion protection	→ See "Approvals"				

Temperature ranges for Ex protection

Temperature class	Ambient temperature
Т6 Т3	$-40 \le Ta \le +40 \text{ °C} [-40 \le Ta \le +104 \text{ °F}]$

Approvals

Approvals included in the scope of delivery

Logo	Description	Country
CE	EU declaration of conformity	European Union
	EMC directive, interference emission (group 1, class B) and immunity per EN 61326-1:2013 (industrial application), EN 61326-2-3:2013 $^{\rm 1)}$	
	Pressure equipment directive	
	RoHS directive	
CHEDC	EHEDG ²⁾ Hygienic Equipment Design	European Community

1) With electrostatic discharge, a short-term, increased error of up to 1 % of the nominal measuring range can occur.

2) EHEDG conformity only in connection with the marked process connections

Optional approvals

Logo	Description	Country	
CE	EU declaration of conformity	European Union	
¢,	ATEX directive Hazardous areas Gas II 2G Ex ia IIC T6 T3 Gb		
	IECEx Hazardous areas Gas Ex ia IIC T6 T3 Gb	International	

Certificates (option)

- 2.2 test report per EN 10204 (e.g. state-of-the-art manufacturing, material proof, indication accuracy)
- 3.1 inspection certificate per EN 10204 (e.g. material proof for wetted metal parts, indication accuracy)
- FDA conformity of the system fill fluid
- 3-A conformity of the diaphragm seal, based on a third party verification
- EHEDG conformity
- Others on request


Patents, property rights

 Diaphragm monitoring for diaphragm seals (US 10794787, NL 2019251)

Approvals and certificates, see website

Dimensions in mm [in]


Clamp connection per DIN 32676, DN 1 1/2"

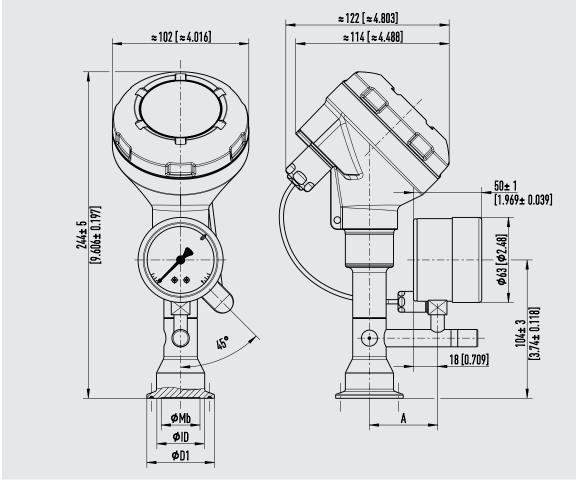

For further information on model UPT-20 see data sheet PE 86.05

14463511.01

Clamp connection per DIN 32676

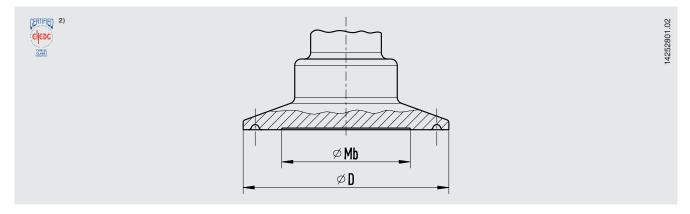

Process connection: Clamp connection per DIN 32676 Pipe standard per DIN 11866 row B and ISO 1127 row 1

DN	PN ¹⁾	Dimensions in mm [in]								
		For pipe outer Ø x wall thickness	Pipe inner Ø	D	Mb					
26.9	40	26.9 x 1.6 [1.059 x 0.063]	23.7 [0.933]	50.5 [1.988]	22 [0.866]					
33.7	40	33.7 x 2 [1.327 x 0.079]	29.7 [1.169]	50.5 [1.988]	25 [0.984]					
42.4	40	42.4 x 2 [1.669 x 0.079]	38.4 [1.512]	64 [2.52]	32 [1.26]					
48.3	40	48.3 x 2 [1.902 x 0.079]	44.3 [1.744]	64 [2.52]	40 [1.575]					
60.3	40	60.3 x 2 [2.374 x 0.079]	56.3 [2.217]	77.5 [3.051]	52 [2.047]					
76.1	25	76.1 x 2 [2.996 x 0.079]	72.1 [2.839]	91 [3.583]	60 [2.362]					

Process connection: Clamp connection per DIN 32676 Pipe standard per DIN 11866 row C or ASME BPE

DN	PN ¹⁾	Dimensions in mm [in]								
		For pipe outer Ø x wall thickness	Pipe inner Ø	D	Mb					
1"	40	25.4 x 1.65 [1 x 0.065]	22.1 [0.87]	50.5 [1.988]	22 [0.866]					
1 1⁄2"	40	38.1 x 1.65 [1.5 x 0.065]	34.8 [1.37]	50.5 [1.988]	32 [1.26]					
2"	40	50.8 x 1.65 [2 x 0.065]	47.5 [1.87]	64 [2.52]	40 [1.575]					
2 ½"	40	63.5 x 1.65 [2.5 x 0.065]	60.2 [2.37]	77.5 [3.051]	52 [2.047]					
3"	25	76.2 x 1.65 [3 x 0.065]	72.9 [2.87]	91 [3.583]	60 [2.362]					


Process connection: Clamp connection per DIN 32676 Pipe standard per BS 4825 part 3 and O.D. tube

DN	PN ¹⁾	Dimensions in mm [in]							
		For pipe outer Ø x wall thickness	Pipe inner Ø	D	Mb				
25.4	40	25.4 x 1.6 [1 x 0.063]	22.2 [0.874]	50.5 [1.988]	22 [0.866]				
38.1	40	38.1 x 1.6 [1.5 x 0.063]	34.9 [1.374]	50.5 [1.988]	32 [1.26]				
50.8	40	50.8 x 1.6 [2 x 0.063]	47.6 [1.874]	64 [2.52]	40 [1.575]				
63.5	40	63.5 x 1.6 [2.5 x 0.063]	60.3 [2.374]	77.5 [3.051]	52 [2.047]				
76.2	25	76.2 x 1.6 [3 x 0.063]	73 [2.874]	91 [3.583]	60 [2.362]				

For maximum pressure range consider pressure rating of clamp.
 EHEDG conformity only in combination with TRI-CLAMP[®] seals Combifit International B.V., The Netherlands.

Further dimensions and higher nominal pressures on request

Clamp connection per DIN 32676

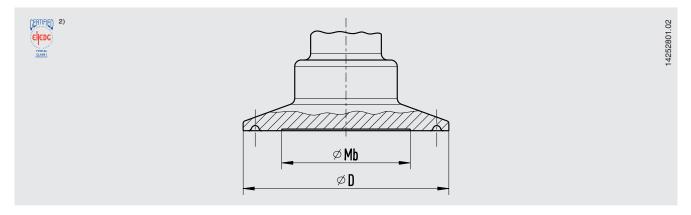


Process connection: Clamp connection per DIN 32676 Pipe standard per DIN 11866 row A and DIN 11850 row 2

DN	PN ¹⁾	Dimensions in mm [in]							
		For pipe outer Ø x wall thickness	Pipe inner Ø	D	Mb				
25	40	29 x 1.5 [x 0.059]	26 [1.024]	50.5 [1.988]	25 [0.984]				
32	40	35 x 1.5 [x 0.059]	32 [1.26]	50.5 [1.988]	29 [1.142]				
40	40	41 x 1.5 [x 0.059]	38 [1.496]	50.5 [1.988]	32 [1.26]				
50	40	53 x 1.5 [x 0.059]	50 [1.969]	64 [2.52]	40 [1.575]				
65	25	70 x 2 [x 0.079]	66 [2.598]	91 [3.583]	59 [2.323]				
80	25	85 x 2 [x 0.079]	81 [3.189]	106 [4.173]	72 [2.835]				
100	25	104 x 2 [x 0.079]	100 [3.937]	119 [4.685]	90 [3.543]				

For maximum pressure range consider pressure rating of clamp.
 EHEDG conformity only in combination with TRI-CLAMP[®] seals Combifit International B.V., The Netherlands.

Further dimensions and higher nominal pressures on request

Clamp connection per ISO 2852

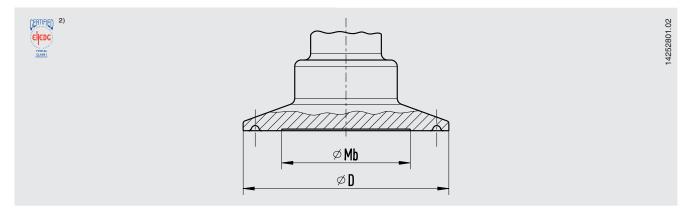

Process connection: Clamp connection per ISO 2852 Pipe standard per ISO 2037 and BS 4825 part 1

DN	PN ¹⁾	Dimensions in mm [in]							
		For pipe outer Ø x wall thickness	er Ø x wall thickness Pipe inner Ø D						
25	40	25 x 1.2 [0.984 x 0.047]	22.6 [0.89]	50.5 [1.988]	22 [0.866]				
33.7	40	33.7 x 1.2 [1.327 x 0.047]	31.3 [1.232]	50.5 [1.988]	25 [0.984]				
38	40	38 x 1.2 [1.496 x 0.047]	35.6 [1.402]	50.5 [1.988]	32 [1.26]				
40	40	40 x 1.2 [1.575 x 0.047]	37.6 [1.48]	64 [2.52]	32 [1.26]				
51	40	51 x 1.2 [2.008 x 0.047]	48.6 [1.913]	64 [2.52]	40 [1.575]				
63.5	40	63.5 x 1.6 [2.5 x 0.063]	60.3 [2.374]	77.5 [3.051]	52 [2.047]				
70	25	70 x 1.6 [2.756 x 0.063]	66.8 [2.63]	91 [3.583]	60 [2.362]				
76.1	25	76.1 x 1.6 [2.996 x 0.063]	72.9 [2.87]	91 [3.583]	60 [2.362]				
88.9	25	88.9 x 2 [3.5 x 0.079]	84.9 [3.343]	106 [4.173]	72 [2.835]				
101.6	25	101.6 x 2 [4 x 0.079]	97.6 [3.843]	119 [4.685]	90 [3.543]				

For maximum pressure range consider pressure rating of clamp.
 EHEDG conformity only in combination with TRI-CLAMP[®] seals Combifit International B.V., The Netherlands.


Further dimensions and higher nominal pressures on request

Aseptic threaded pipe connection per DIN 11864-1



With threaded coupling

With liner and union nut

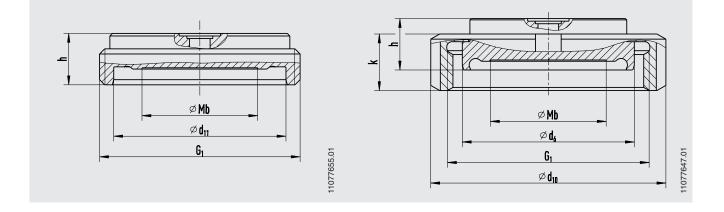

Type of process connection: Aseptic threaded pipe connection per DIN 11864-1 form A Pipe standard per DIN 11866 row A or DIN 11850 row 2


DN	For pipe	PN ¹⁾	Dimensi	ons in mn	ո [in]					
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	G ₁	h	d ₁₀	k	Aseptic O-ring
25	29 x 1.5 [1.142 x 0.071]	40	22 [0.866]	42.9 [1.689]	43 [1.693]	RD 52 x ½ [2.047 x ½]	20 [0.787]	63 [2.48]	21 [0.827]	28 x 3.5 [1.102 x 0.1378]
32	35 x 1.5 [1.378 x 0.071]	40	25 [0.984]	48.9 [1.925]	49 [1.929]	RD 58 x ½ [2.283 x ½]	20 [0.787]	70 [2.756]	21 [0.827]	34 x 5 [1.339 x 0.197]
40	41 x 1.5 [1.614 x 0.071]	40	35 [1.378]	54.9 [2.161]	55 [2.165]	RD 65 x ½ [2.559 x ½]	20 [0.787]	78 [3.071]	21 [0.827]	40 x 5 [1.575 x 0.197]
50	53 x 1.5 [2.087 x 0.071]	25	45 [1.772]	66.9 [2.634]	67 [2.638]	RD 78 x ½ [3.071 x ½]	20 [0.787]	92 [3.622]	22 [0.866]	52 x 5 [2.047 x 0.197]
65	70 x 2 [2.756 x 0.078]	25	60 [2.362]	84.9 [3.343]	85 [3.346]	RD 95 x ½ [3.74 x ½]	20 [0.787]	112 [4.409]	25 [0.984]	68 x 5 [2.677 x 0.197]
80	85 x 2 [3.346 x 0.078]	25	72 [2.835]	98.9 [3.894]	99 [3.898]	RD 110 x ¼ [4.331 x ¼]	20 [0.787]	127 [5]	29 [1.142]	83 x 5 [3.268 x 0.197]
100	104 x 2 [4.094 x 0.078]	25	90 [3.543]	118.9 [4.681]	119 [4.685]	RD 130 x ¼ [5.118 x ¼]	20 [0.787]	148 [5.827]	31 [1.22]	102 x 5 [4.016 x 0.197]

Aseptic threaded pipe connection per DIN 11864-1



With threaded coupling

With liner and union nut

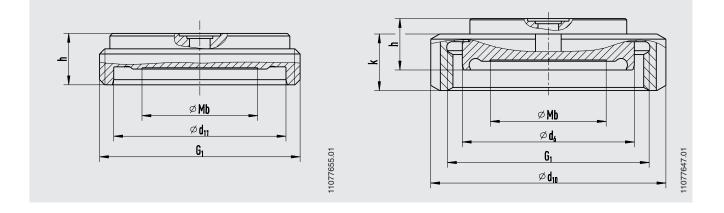

Type of process connection: Aseptic threaded pipe connection per DIN 11864-1 form A Pipe standard per DIN 11866 row B or DIN ISO 1127 row 1


DN	For pipe	PN 1)	Dimensi	ons in m	m [in]					
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	G ₁	h	d ₁₀	k	Aseptic O-ring
26.9	26.9 x 1.6 [1.059 x 0.063]	40	22 [0.866]	42.9 [1.689]	43 [1.693]	RD 52 x ½ [2.047 x ½]	20 [0.787]	63 [2.48]	21 [0.827]	26 x 3.5 [1.024 x 0.1378]
33.7	33.7 x 2 [1.327 x 0.078]	40	25 [0.984]	48.9 [1.925]	49 [1.929]	RD 58 x ½ [2.283 x ½]	20 [0.787]	70 [2.756]	21 [0.827]	32 x 5 [1.26 x 0.197]
42.4	42.4 x 2 [1.669 x 0.078]	25	35 [1.378]	54.9 [2.161]	55 [2.165]	RD 65 x ½ [2.559 x ½]	20 [0.787]	78 [3.071]	21 [0.827]	40.5 x 5 [1.594 x 0.197]
48.3	48.3 x 2 [1.902 x 0.078]	25	45 [1.772]	66.9 [2.634]	67 [2.638]	RD 78 x ½ [3.071 x ½]	20 [0.787]	92 [3.622]	22 [0.866]	46.5 x 5 [1.831 x 0.197]
60.3	60.3 x 2 [2.374 x 0.078]	25	60 [2.362]	84.9 [3.343]	85 [3.346]	RD 95 x ¼ [3.74 x ¼]	20 [0.787]	112 [4.409]	25 [0.984]	58.5 x 5 [1.831 x 0.197]
76.1	76.1 x 2 [2.996 x 0.078]	25	72 [2.835]	98.9 [3.894]	99 [3.898]	RD 110 x ¼ [4.331 x ¼]	20 [0.787]	127 [5]	29 [1.142]	73.5 x 5 [2.894 x 0.197]
88.9	88.9 x 2.3 [3.5 x 0.091]	25	90 [3.543]	118.9 [4.681]	119 [4.685]	RD 130 x ¼ [5.118 x ¼]	20 [0.787]	148 [5.827]	31 [1.22]	86.5 x 5 [3.406 x 0.197]

Aseptic threaded pipe connection per DIN 11864-1



With threaded coupling

With liner and union nut

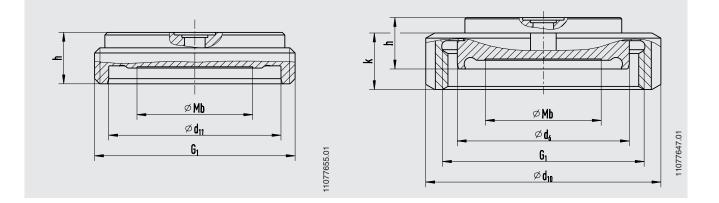

Type of process connection: Aseptic threaded pipe connection per DIN 11864-1 form A Pipe standard per DIN 11866 row C or ASME BPE 1997


DN	For pipe	PN 1)	Dimensi	mensions in mm [in]						
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d11	G1	h	d ₁₀	k	Aseptic O-ring
1"	25.4 x 1.65 [1.831 x 0.065]	40	22 [0.866]	42.9 [1.689]	43 [1.693]	RD 52 x ½ [2.047 x ½]	20 [0.787]	63 [2.48]	21 [0.827]	24 x 3,5 [0.945 x 0.1378]
1 1⁄2"	42.4 x 1.65 [1.669 x 0.065]	40	32 [1.26]	54.9 [2.161]	55 [2.165]	RD 65 x ½ [2.559 x ½]	20 [0.787]	78 [3.071]	21 [0.827]	37 x 5 [1.457 x 0.197]
2"	48.3 x 1.65 [1.902 x 0.065]	25	45 [1.772]	66.9 [2.634]	67 [2.638]	RD 78 x ½ [3.071 x ½]	20 [0.787]	92 [3.622]	22 [0.866]	50 x 5 [1.969 x 0.197]
2 ½"	60.3 x 1.65 [2.374 x 0.065]	25	52 [2.047]	84.9 [3.343]	85 [3.346]	RD 95 x ¼ [3.74 x ½]	20 [0.787]	112 [4.409]	25 [0.984]	62 x 5 [2.441 x 0.197]
3"	76.1 x 1.65 [2.996 x 0.065]	25	60 [2.362]	98.9 [3.894]	99 [3.898]	RD 110 x ¼ [4.331 x ¼]	20 [0.787]	127 [5]	29 [1.142]	75 x 5 [2.953 x 0.197]
4"	88.9 x 2.11 [3.5 x 0.083]	25	90 [3.543]	118.9 [4.681]	119 [4.685]	RD 130 x ¼ [5.118 x ¼]	20 [0.787]	148 [5.827]	31 [1.22]	100 x 5 [3.937 x 0.197]

Aseptic flange connection per DIN 11864-2



With flange with groove

With flange with notch

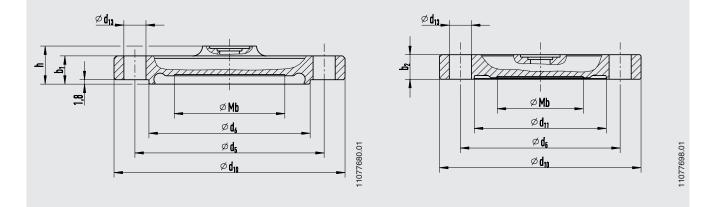

Type of process connection: Aseptic flange connection per DIN 11864-2 form A Pipe standard per DIN 11866 row A or DIN 11850 row 2


DN	For pipe	PN ¹⁾	Dimens	sions in	mm [in]							
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	d ₅	d ₁₀	h	b ₁	b ₂	d ₁₃	Aseptic O-ring
25	29 x 1.5 [1.142 x 0.071]	25	22 [0.866]	38.3 [1.508]	38.4 [1.512]	53 [2.087]	70 [2.756]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	28 x 3.5 [1.102 x 0.1378]
32	35 x 1.5 [1.378 x 0.071]	25	25 [0.984]	47.6 [1.878]	47.7 [1.878]	59 [2.323]	76 [2.992]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	34 x 5 [1.339 x 0.197]
40	41 x 1.5 [1.614 x 0.071]	25	35 [1.378]	53.6 [2.11]	53.7 [2.114]	65 [2.559]	82 [3.228]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	40 x 5 [1.575 x 0.197]
50	53 x 1.5 [2.087 x 0.071]	16	45 [1.772]	65.6 [2.583]	65.7 [2.587]	77 [3.032]	94 [3.7]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	52 x 5 [2.047 x 0.197]
65	70 x 2 [2.756 x 0.078]	16	60 [2.362]	81.6 [3.213]	81.7 [3.217]	95 [3.74]	113 [4.449]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	8 x Ø 9 [0.354]	68 x 5 [2.677 x 0.197]
80	85 x 2 [3.346 x 0.078]	16	72 [2.835]	97.6 [3.843]	97.7 [3.846]	112 [4.409]	133 [5.236]	17.5 [0.689]	13.5 [0.531]	12 [0.472]	8 x Ø 11 [0.433]	83 x 5 [3.268 x 0.197]
100	104 x 2 [4.094 x 0.078]	16	90 [3.543]	116.6 [4.591]	116.7 [4.594]	137 [5.394]	159 [6.2598]	19.5 [0.768]	15.5 [0.61]	14 [0.551]	8 x Ø 11 [0.433]	102 x 5 [4.016 x 0.197]

Aseptic flange connection per DIN 11864-2



With flange with groove

With flange with notch

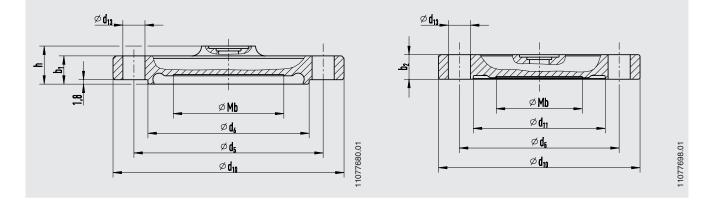

Type of process connection: Aseptic flange connection per DIN 11864-2 form A Pipe standard per DIN 11866 row B or DIN ISO 1127 row 1


DN	For pipe	PN ¹⁾	Dimens	ions in m	ım [in]							
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	d ₅	d ₁₀	h	b ₁	b ₂	d ₁₃	Aseptic O-ring
26.9	26.9 x 1.6 [1.059 x 0.063]	25	22 [0.866]	36 [1.417]	36.1 [1.421]	52 [2.047]	69 [2.717]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	26 x 3.5 [1.024 x 0.1378]
33.7	33.7 x 2 [1.327 x 0.078]	25	25 [0.984]	45.3 [1.783]	45.4 [1.787]	57 [2.244]	74 [2.913]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	32 x 5 [1.2598 x 0.197]
42.4	42.4 x 2 [1.669 x 0.078]	16	35 [1.378]	54 [2.126]	54.1 [2.1299]	65 [2.559]	82 [3.228]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	40.5 x 5 [1.594 x 0.197]
48.3	48.3 x 2 [1.902 x 0.078]	16	40 [1.575]	59.9 [2.358]	60 [2.362]	71 [2.795]	88 [3.465]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	46.5 x 5 [1.831 x 0.197]
60.3	60.3 x 2 [2.374 x 0.078]	16	52 [2.047]	71.9 [2.831]	72 [2.835]	85 [3.346]	103 [4.055]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	8 x Ø 9 [0.354]	58.5 x 5 [1.831 x 0.197]
76.1	76.1 x 2 [2.996 x 0.078]	16	60 [2.362]	88.1 [3.469]	88.1 [3.469]	104 [4.094]	125 [4.921]	17.5 [0.689]	13.5 [0.531]	12 [0.472]	8 x Ø 11 [0.433]	73.5 x 5 [2.894 x 0.197]
88.9	88.9 x 2.3 [3.5 x 0.091]	16	72 [2.835]	100.9 [3.972]	101 [3.976]	116 [4.567]	137 [5.394]	17.5 [0.689]	13.5 [0.531]	12 [0.472]	8 x Ø 11 [0.433]	86.5 x 5 [3.406 x 0.197]

Aseptic flange connection per DIN 11864-2



With flange with groove

With flange with notch

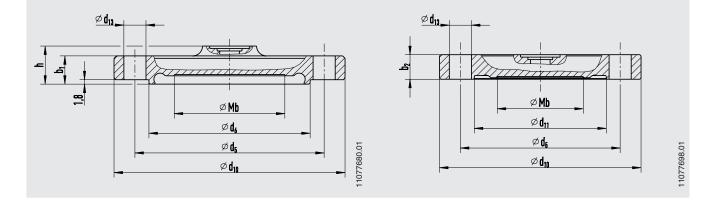

Type of process connection: Aseptic flange connection per DIN 11864-2 form A Pipe standard per DIN 11866 row C or ASME BPE 1997


DN	For pipe	PN 1)	Dimens	ions in	mm [in]							
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	G ₁	d ₁₀	h	b ₁	b ₂	d ₁₃	Aseptic O-ring
1"	25.4 x 1.65 [1.831 x 0.065]	25	22 [0.866]	34.3 [1.354]	34.4 [1.354]	49 [1.929]	66 [2.598]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	24 x 3.5 [0.945 x 0.1378]
1 1⁄2"	42.4 x 1.65 [1.669 x 0.065]	25	32 [1.26]	50.4 [1.984]	50.4 [1.984]	62 [2.44]	79 [3.11]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	37 x 5 [1.457 x 0.197]
2"	48.3 x 1.65 [1.902 x 0.065]	16	45 [1.772]	63 [2.48]	63 [2.48]	75 [2.953]	92 [3.622]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	50 x 5 [1.969 x 0.197]
2 ½"	60.3 x 1.65 [2.374 x 0.065]	16	52 [2.047]	75.8 [2.984]	75.9 [2.988]	89 [3.504]	107 [4.213]	15.5 [0.61]	11.5 [0.071]	10 [0.394]	4 x Ø 9 [0.354]	62 x 5 [2.441 x 0.197]
3"	76.1 x 1.65 [2.996 x 0.065]	16	60 [2.362]	89.5 [3.524]	89.6 [3.528]	104 [4.094]	125 [4.921]	17.5 [0.689]	13.5 [0.531]	12 [0.472]	8 x Ø 11 [0.433]	75 x 5 [2.953 x 0.197]
4"	88.9 x 2.11 [3.5 x 0.083]	16	90 [3.543]	114.2 [4.496]	114.3 [4.5]	135 [5.315]	157 [6.181]	19.5 [0.768]	15.5 [0.61]	14 [0.551]	8 x Ø 11 [0.433]	100 x 5 [3.937 x 0.197]

Aseptic clamp connection per DIN 11864-3



With clamp with groove

With clamp with notch

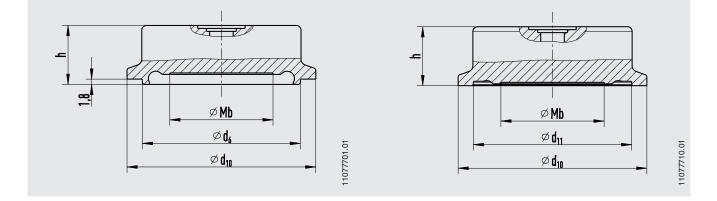

Type of process connection: Aseptic clamp connection DIN 11864-3 form A Pipe standard per DIN 11866 row A or DIN 11850 row 2


DN	For pipe	PN 1)	Dimensions i	n mm [in]				
	Outer Ø x wall thick- ness in mm [in]		Mb	d ₆	d ₁₁	h	d ₁₀	Aseptic O-ring
25	29 x 1.5 [1.142 x 0.071]	40	22 [0.866]	38.3 [1.508]	38.4 [1.512]	20 [0.787]	50.5 [1.988]	28 x 3.5 [1.102 x 0.1378]
32	35 x 1.5 [1.378 x 0.071]	40	25 [0.984]	47.6 [1.878]	47.7 [1.878]	20 [0.787]	50.5 [1.988]	34 x 5 [1.339 x 0.197]
40	41 x 1.5 [1.614 x 0.071]	40	35 [1.378]	53.6 [2.11]	53.7 [2.114]	20 [0.787]	64 [2.5197]	40 x 5 [1.575 x 0.197]
50	53 x 1.5 [2.087 x 0.071]	25	45 [1.772]	65.6 [2.583]	65.7 [2.587]	20 [0.787]	77.5 [3.051]	52 x 5 [2.047 x 0.197]
65	70 x 2 [2.756 x 0.078]	25	60 [2.362]	81.6 [3.213]	81.7 [3.217]	20 [0.787]	91 [3.583]	68 x 5 [2.677 x 0.197]
80	85 x 2 [3.346 x 0.078]	16	72 [2.835]	97.6 [3.843]	97.7 [3.846]	20 [0.787]	106 [4.173]	83 x 5 [3.268 x 0.197]
100	104 x 2 [4.094 x 0.078]	16	90 [3.543]	116.6 [4.591]	116.7 [4.594]	20 [0.787]	130 [5.118]	102 x 5 [4.016 x 0.197]

Aseptic clamp connection per DIN 11864-3



With clamp with groove

With clamp with notch

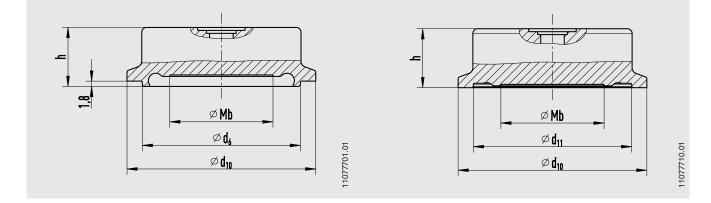

Type of process connection: Aseptic clamp connection DIN 11864-3 form A Pipe standard per DIN 11866 row B or DIN ISO 1127 row 1


DN	For pipe	PN ¹⁾	Dimensions	in mm [in]				
	Outer Ø x wall thick- ness in mm [in]		Мb	d ₆	d ₁₁	h	d ₁₀	Aseptic O-ring
26.9	26.9 x 1.6 [1.059 x 0.063]	40	22 [0.866]	36.0 [1.417]	36.1 [1.421]	20 [0.787]	50.5 [1.988]	26 x 3.5 [1.024 x 0.1378]
33.7	33.7 x 2 [1.327 x 0.078]	40	25 [0.984]	45.3 [1.783]	45.3 [1.783]	20 [0.787]	50.5 [1.988]	32 x 5 [1.2598 x 0.197]
42.4	42.4 x 2 [1.669 x 0.078]	25	35 [1.378]	54.0 [2.126]	54.1 [2.1299]	20 [0.787]	64 [2.5197]	40.5 x 5 [1.594 x 0.197]
48.3	48.3 x 2 [1.902 x 0.078]	25	40 [1.575]	59.9 [2.358]	60 [2.362]	20 [0.787]	64 [2.5197]	46.5 x 5 [1.831 x 0.197]
60.3	60.3 x 2 [2.374 x 0.078]	25	52 [2.047]	71.9 [2.831]	72.0 [2.835]	20 [0.787]	91 [3.583]	58.5 x 5 [2.303 x 0.197]
76.1	76.1 x 2 [2.996 x 0.078]	16	60 [2.362]	88.1 [3.469]	88.2 [3.472]	20 [0.787]	106 [4.173]	73.5 x 5 [2.894 x 0.197]
88.9	88.9 x 2 [3.5 x 0.078]	16	72 [2.835]	100.9 [3.972]	101 [3.976]	25 [0.984]	119 [4.685]	86.5 x 5 [3.406 x 0.197]

Aseptic clamp connection per DIN 11864-3



With clamp with groove


With clamp with notch

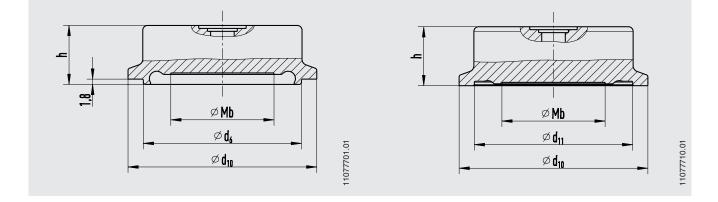

Type of process connection: Aseptic clamp connection DIN 11864-3 form A Pipe standard per DIN 11866 row C or ASME BPE 1997

DN	For pipe	PN 1)	Dimensions i	n mm [in]				
	Outer Ø x wall thickness in mm [in]		Mb	d ₆	d ₁₁	h	d ₁₀	Aseptic O-ring
1"	25.4 x 1.65 [1.831 x 0.065]	40	22 [0.866]	34.3 [1.35]	34.4 [1.354]	20 [0.787]	50.5 [1.988]	24 x 3.5 [0.945 x 0.1378]
1 1⁄2"	42.4 x 1.65 [1.669 x 0.065]	40	32 [1.26]	50.4 [1.984]	50.5 [1.988]	20 [0.787]	64 [2.5197]	37 x 5 [1.457 x 0.197]
2"	48.3 x 1.65 [1.902 x 0.065]	25	45 [1.772]	63 [2.48]	63 [2.48]	20 [0.787]	77.5 [3.051]	50 x 5 [1.969 x 0.197]
2 ½"	60.3 x 1.65 [2.374 x 0.065]	25	52 [2.047]	75.8 [2.984]	75.9 [2.988]	20 [0.787]	91 [3.583]	62 x 5 [2.441 x 0.197]
3"	76.1 x 1.65 [2.996 x 0.065]	16	60 [2.362]	89.5 [3.524]	89.6 [3.528]	20 [0.787]	106 [4.173]	75 x 5 [2.953 x 0.197]
4"	88.9 x 2.11 [3.5 x 0.083]	16	90 [3.543]	114.2 [4.496]	114.3 [4.5]	25 [0.984]	130 [5.118]	100 x 5 [3.937 x 0.197]


Ingold connection with union nut

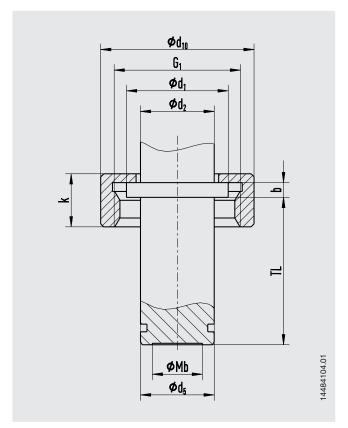

DN	N PN Dimensions in mm [in]									
		Mb	d 5	d ₁	d ₂	b	TL	G ₁	d ₁₀	k
25	25	17 [0.669]	25 [0.984]	34.5 [1.358]	25 [0.984]	5 [0.917]	50 [1.969]	G 1 ¼	52 [2.047]	18 [0.709]
40	25	29 [1.142]	40 [1.575]	55 [2.165]	25 [0.984]	5 [0.917]	56 [2.205]	G 2	78 [3.071]	27 [1.063]
50	25	38 [1.496]	50 [1.969]	55.5 [2.185]	25 [0.984]	5 [0.917]	60 [2.362]	G 2	78 [3.071]	32 [1.26]

This process connection is delivered with a sealing from NBR, FKM or EPDM.

Diaphragm monitoring via DTM

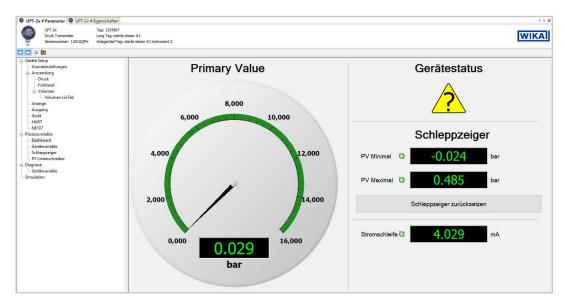

For HART[®] output signals, a DTM is available in accordance with the FDT standard. The DTM provides a self-explanatory and clear user interface for all setup and control processes of the transmitter. For testing purposes, it is also possible to simulate process values and archive the parameter data.

Recording of the measured values is available for diagnostic purposes.


Diagnostics: Diaphragm rupture

In the event of a diaphragm rupture in the first diaphragm, the instrument status will be set to "Diaphragm rupture". This allows the user to analyse faults to the second, while the instrument continues to monitor the process pressure without restriction. The user thus has the advantage of reducing to a minimum any production errors that may occur. No contamination of the process medium or the environment has occurred at this time.



Tank volume calculation

The additional tank volume calculation of the DTM function can be used to reproduce any optional tank geometry. The corresponding linearisation table is generated automatically. The linearisation table can be transferred directly to the transmitter.

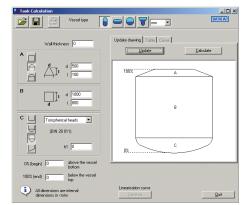

Accessories and spare parts

Description		Order number		
i i i i i i i i i i i i i i i i i i i	HART® modem for USB interface, specifically designed for use with notebooks (model 010031)	11025166		
	HART [®] modem for RS-232 interface (model 010001)	7957522		
March	HART® modem for Bluetooth® interface Ex ia IIC (model 010041)	11364254		
V 🕈 🖓	PowerXpress HART [®] modem, with optional power supply (model 010031P) 1			
	Overvoltage protection for transmitters, 4 20 mA, M20 x 1.5, series connection	14002489		
	Display and operating unit, model DI-PT-U The display and operating unit can be attached in 90° steps. The display and operating unit features a main display and an additional display. The main display shows the output signal. The additional display shows different values, at the same time as the main display - these values can be selected by the user. The process pressure transmitter can be configured through the display and operating unit. Only this display may be used for installation into the process transmitter.	14090181		
	Hygienic cable gland M20 x 1.5 Cable diameter: 6 12 mm [0.24 0.47 in]	11348691		

Instruments for on-site calibration

Model	Description
СРС-КІ	 Pneumatic service kit, accuracy 0.1 % FS (0.05 % FS or 0.025 % FS also available) Model CPG1500 precision digital pressure gauge Model CPP30 pneumatic hand test pump, pressure generation -0.95 +35 bar Adapter set Service case → see data sheet CT 93.01
СРН700	 Portable process calibrator, accuracy 0.025 % FS Model CPH7000 process calibrator, integrated manual pressure generation -0.85 +25 bar Power supply unit Service case → see data sheet CT 15.51
CPH76	 Portable pressure calibrator, accuracy 0.025 % FS Model CPH7650 pressure calibrator, integrated electrical pressure generation -0.85 +20 bar Test cables Battery charger → see data sheet CT 17.02

FS = Full span = end of measuring range - start of measuring range

Calibration adapter

Description	Order number
Calibration adapter TRI-CLAMP®, 1 1/2"	11563206
Calibration adapter TRI-CLAMP®, 2"	14332415

Other calibration adapters on request

WIKA-Cal calibration software

Easy and fast creation of a high-quality calibration certificate

The WIKA-Cal calibration software is used for generating calibration certificates or logger protocols for pressure measuring instruments and is available as a demo version for a cost-free download.

A template helps the user and guides him through the creation process of a document.

In order to switch from the demo version to a full version of the respective template, a USB stick with the template must be purchased.

The pre-installed demo version automatically changes to the selected full version when the USB stick is inserted and remains available as long as the USB stick is connected to the computer.

- Creation of calibration certificates for mechanical and electronic pressure measuring instruments
- A calibration assistant guides you through the calibration
- Automatic generation of the calibration steps
- Generation of 3.1 certificates per DIN EN 10204
- Creation of logger protocols
- User-friendly interface
- Languages: German, English, Italian and more due with software updates

For further information see data sheet CT 95.10



Calibration certificates can be created with the Cal-Template and logger protocols can be created with the Log-Template.

Cal
Demo

Cal Demo

Generation of calibration certificates limited to 2 measuring points, with automatic initiation of pressures via a pressure controller.



Cal Light

Generation of calibration certificates with no limitations on measuring points, without automatic initiation of pressures via a pressure controller.



Log Demo

Creation of data logger test reports, limited to 5 measured values.



Log

Creation of data logger test reports without limiting the measured values.

Calibration certi Kaliniamoph	fcate		WIKA
		Cartilizata No. Jaupris No.	1000-20
Carenal Nijeran Santara Santar	No alian Kasau Kasau Kasau Kasau	Denins under Verd (Antreapperstein) für Reinersteiner Rei	NT) N An An An An An An An An An An An An An
Endersonantial constit impetion production of		Para disalitation internationag	
a preserve	100.001		110000-000-00
· material	100.00		
. 513555	181.175		
Newsrament scralit Newsramperper	-		
The local division in which the local division in which the local division in the local			
-	Auto Die		
Table State State	Salar Ville		
Same areas	10 - 10 Mar 1		
The second second			



Ordering information

Model / Case head / Surface roughness / System fill fluid / Measuring range / Accuracy / Output signal / Electrical connection / Process connection / Certificates

© 05/2021 WIKA Alexander Wiegand SE & Co. KG, all rights reserved. The specifications given in this document represent the state of engineering at the time of publishing. We reserve the right to make modifications to the specifications and materials.

WIKA data sheet DS 95.11 · 05/2021



Your WIKA Sales Partner

ICS Schneider Messtechnik GmbH Briesestrasse 59 D-16562 Hohen Neuendorf / OT Bergfelde Tel.: +49 3303 5040-66 Fax: +49 3303 5040-68 E-Mail: info@ics-schneider.de



WIKA Alexander Wiegand SE & Co. KG Alexander-Wiegand-Straße 30 63911 Klingenberg/Germany Tel. +49 9372 132-0 Fax +49 9372 132-406 info@wika.de www.wika.de

Page 27 of 27